

Mix it well, for best results.

DILUTION CHART FOR 5 LITRE CONTAINERS WITH PUMP



SIZE OF CONTAINER TO BE FILLED

DILUTION	500 ml	750 ml	1 Litre	5 Litre	10 Litre	20 Litre
1:10 / 10% Discard jars, TC & body spills	5 pumps or 50 ml	Self-dose or 75 ml	10 pumps or 100 ml	50 pumps or 500 ml	-	-
1:50 / 2% High risk disinfection of organisms in dirty conditions	1 pump or 10 ml	Self-dose or 15 ml	2 pumps or 20 ml	5 pumps or 100 ml	20 pumps or 200 ml	40 pumps or 400 ml
1:100 / 1% General, intermediate risk use	Self-dose or 5 ml	Self-dose or 7.5 ml	1 pump or 10 ml	5 pumps or 50 ml	10 pumps or 100 ml	20 pumps or 200 ml



Chemgene MEDLAB Multi-Surface Disinfectant

Clear, Unfragranced Art. No. **SKU075A**

Blue, Eucalyptus Art. No. **SKU077A**

Pump, 10 ml dose Art. No. **TR021**

FIND OUT MORE AT

 starlab.click/disinfectant