This is your **Microplate Mixer** user manual.

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1	Safety instructions	4
2	Product description	4
3	Inspection	4
4	Ambient conditions	5
5	Operation	5
5.1	Control elements	5
5.2	Control	5
5.3	Speed range	5
6	Maintenance and cleaning	6
6.1	Warranty	6
6.2	Service	6
7	Technical specifications	7
8	Ordering information	7
8.1	Accessories	7

Latest version of the Declaration of Conformity is available at:
starlab.click/doc-microplatemixer

1 Safety instructions

Before operation

- › Read the instruction manual carefully before use.
- › Ensure that only trained staff use the instrument.
- › Before each use, check both the instrument and the accessories for damage. Do not use damaged components.
- › Make sure the socket is earthed before use.
- › The operation voltage noted on the label must match.
- › Set the instrument on a stable, clean, dry, non-slip and fireproof surface.
- › Do not operate in hazardous areas.
- › Accessories must be securely attached.
- › Always disconnect the plug before fitting accessories.
- › Safe operation is only guaranteed with the original accessories from Starlab. You will find the complete list in chapter 8.
- › When working, wear personal protection to avoid the risk from splashing of liquids and possible glass breakage which can result from mechanical vibrations.
- › The instrument must only be opened by a qualified and trained technician.

During operation

- › Set the "speed-control knob" to the lowest position first and then increase the speed.
- › Reduce the speed if medium splashes out of the vessel or the instrument is unstable on the bench.
- › Keep a minimum distance of 100 mm clear space around the instrument and away from walls and other instruments.

2 Product description

The instrument is designed for mixing liquids.

Normal use: The instrument is designed for use in professional laboratories. Do not use it in residential areas.

3 Inspection

Unpack the equipment carefully and check for any damages. If the instrument is faulty you should contact Starlab or your Starlab distributor promptly.

The packing includes:

- › One Microplate Mixer N2400-8040
- › One EU-power supply and one UK-power supply
- › One instruction manual

Please check all items are present.

4 Ambient conditions

- › Power supply: 100 – 240 V 50/60Hz
- › Ambient temperature: 4 – 40 °C
- › Relative humidity: < 80 %

Place the instrument on a safe and stable surface.

5 Operation

5.1 Control elements

5.2 Control

- › Make sure that the required operating voltage and power supply voltage match.
- › Plug in the power adapter.
- › Switch on the device at the power switch.
- › Set the rated mix speed and the device begins to work.
- › Switch off the device at the power switch.

5.3 Speed range

- › Attachment for one microtiter plate: 0 – 1500 rpm
- › Attachment for two microtiter plates: 0 – 1000 rpm
- › Tube adapters: 0 – 1500 rpm

6 Maintenance and cleaning

- › Proper maintenance and cleaning will keep the instrument in good working order and will help extend its life time.
- › Wear protective gloves during cleaning of the instrument.
- › Be careful not spray cleaning fluid into the instrument. Unplug the power cord when cleaning.
- › Please use only Isopropyl alcohol and water containing tenside. Before using other methods or cleaning agents the user must ascertain with the manufacturer that this will not destroy the instrument.
- › When not in use for long periods, switch off the instrument and store it in a dry, clean and stable place at room temperature.

6.1 Warranty

This instrument is under warranty to be free from defects in materials and workmanship under normal use and service, for a period of 24 months from the date of invoice. The warranty claim expires if the instrument is opened by untrained or unqualified staff. Our “General terms and conditions” apply.

6.2 Service

If help is needed, please contact Starlab or your Starlab distributor. See www.starlabgroup.com for your local contact. Please provide the customer care representative with the following information:

- › Serial Number (on the rear panel)
- › Description of problem (i.e., hardware or software)
- › Methods and procedures adopted to resolve the problems
- › Your contact information

7 Technical specifications

Parameter	Data
Voltage	100 – 240 V
Frequency	50/60 Hz
Power	20 W
Mixing movement	Orbital
Orbital diameter	4.5 mm
Max. mixing weight (with attachments)	0.5 kg
Motor type	Brushless motor
Motor rating input	18 W
Motor rating output	10 W
Permissible ON time	100 %
Speed range	0 – 1500 rpm
Speed display	Scale
Dimensions (WxDxH)	150 x 260 x 80 mm
Weight	3 kg
Permissible ambient temperature	4 – 40 °C
Permissible relative humidity	< 80 %
Protection class acc. to DIN 60529	IP21

8 Ordering information

Article number	Description
N2400-8040	Microplate Mixer

8.1 Accessories

Article number	Description
N2400-8041	Replacement attachment for one microtiter plate
N2400-8042	Attachment for two microtiter plates
N2400-8043	Microplate Mixer Adapter for use with N2400-6012 & N2400-6013
N2400-6012	Tube adapter for 15 x 1.5/2.0 ml Tubes
N2400-6013	Tube adapter for 8 x 15 ml Centrifuge Tubes

Please visit our website www.starlabgroup.com, for details on our other instruments and our comprehensive range of consumables for the laboratory, including PlateOne® deepwell and microplates.

User manual	3
Bedienungsanleitung	9
Instruction initiale	15
Istruzioni d'uso	21

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